

Notes on Using PFV 4.4.0.0 in a Norton Environment


There is a possibility that Norton may detect PFV and, due to an issue on Norton's side, trigger a **BSOD (Blue Screen of Death)**.

This document explains how to exclude PFV from Norton scans to help prevent such incidents.

Step 1: Launch Norton and Click "Security"

From the menu on the left, select **"Security"**.

This section allows you to manage virus scans and exclusion settings.

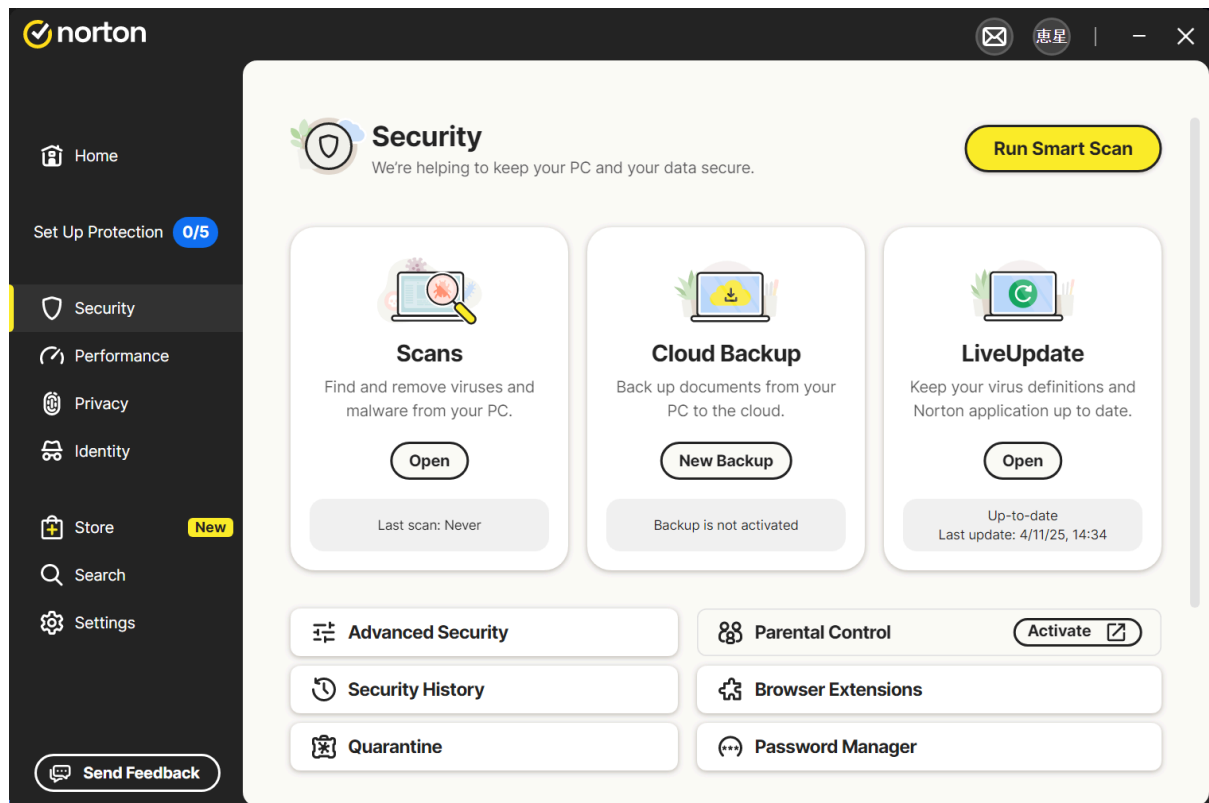


Step 2: Click "Scans"

Under the "Security" tab, click the **"Open"** button under **"Scans"**.

Step 3: Open the "Exclusions" Tab

From the top tab menu, select **"Exclusions"**.

This screen allows you to specify files or folders to exclude from scans.

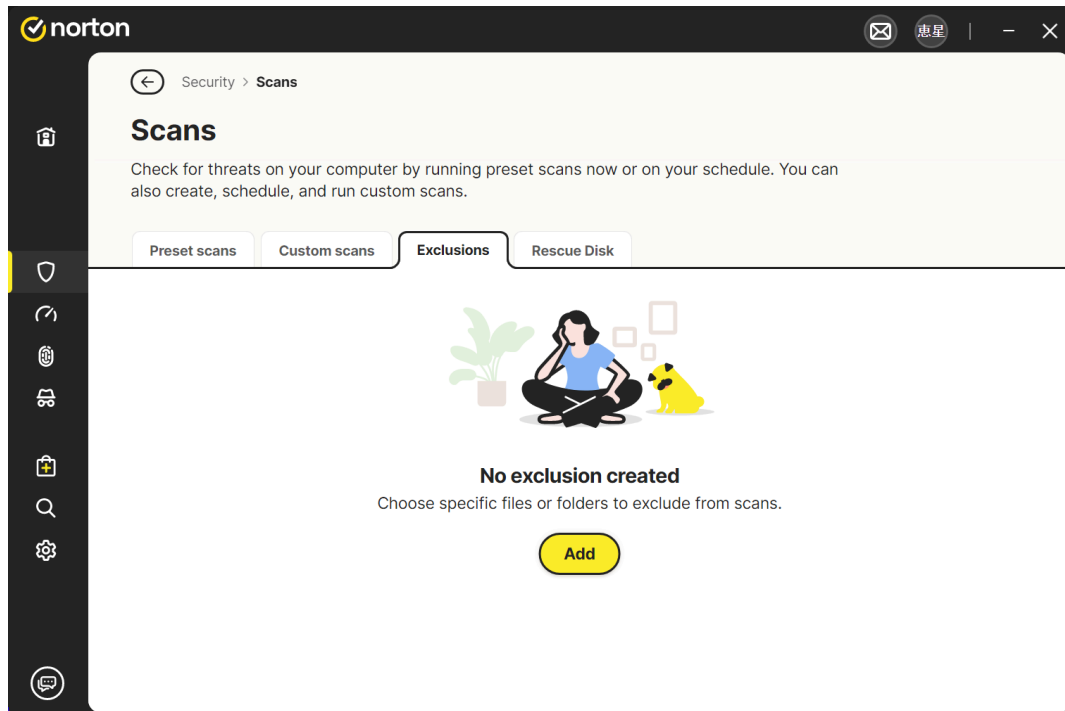

Step 4: Click "Add" and Select the Folder to Exclude


Click the **"Add"** button and select the following folder:

Unset

`C:\Program Files\Photron\Photron FASTCAM Viewer 4\`

Step 5: Confirm the Folder Has Been Added

Make sure the selected folder appears in the list.
If you see it, the setup is complete

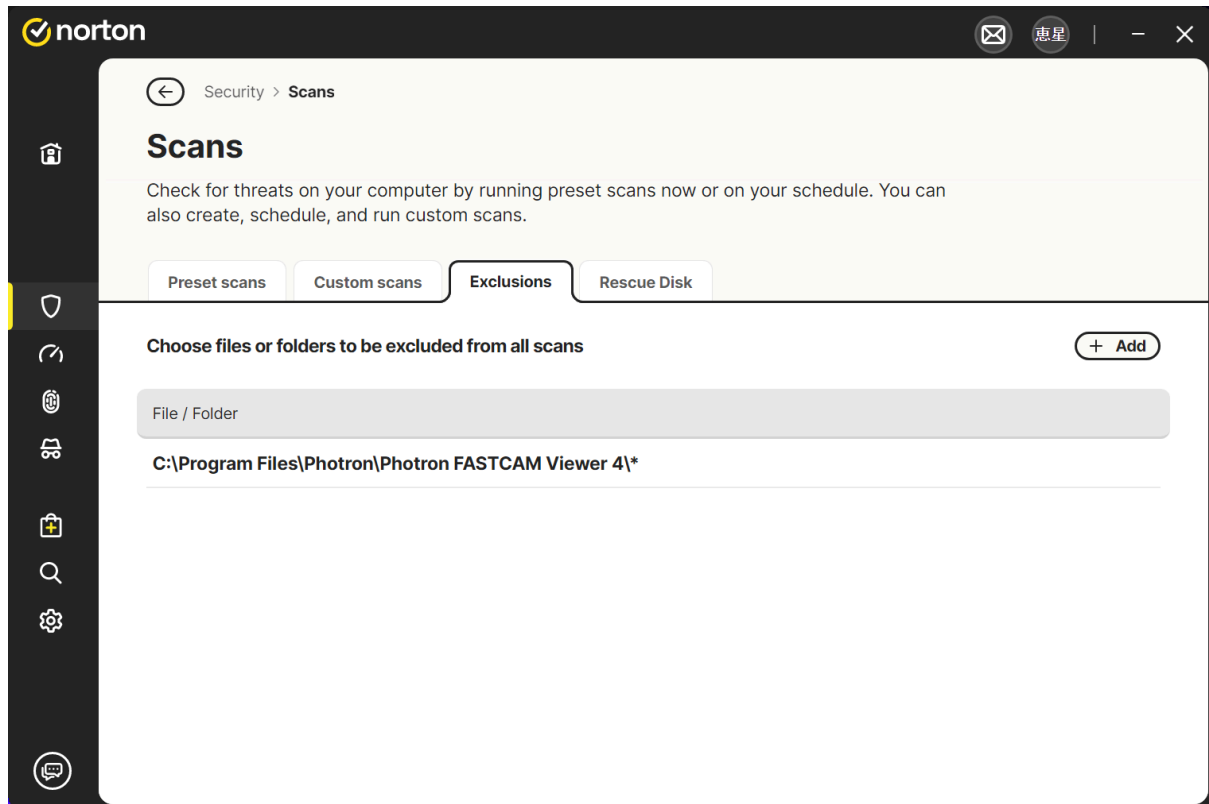

Step 6: Scan the PFV4 Folder

Open **File Explorer** and navigate to the following path:

`C:\Program Files\Photron\Photron FASTCAM Viewer 4\`

Right-click the folder, select **Norton**, then left-click "**Scan Now**" or "**Scan Selected Items**".